# Levenhuk MED 5M Microscope Digital Camera with 9.4" LCD Screen

**EN** User manual

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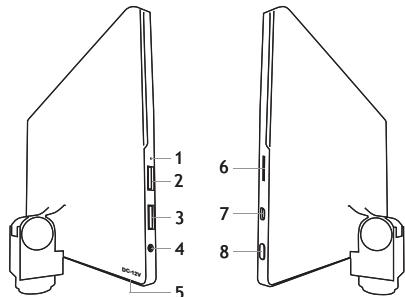
Never look at the sources of bright light or lasers through the camera or the microscope, as this may cause permanent eye damage and even blindness.

## General information

Levenhuk MED Digital Camera is a professional accessory for trinocular microscopes. Its 9.4" color LCD touch screen and 5Mpx digital camera allow you to conduct high-precision research in real time, and save images and video for further use.

### Parts of digital camera:

1. Microphone
2. USB port
3. USB mouse port
4. Earphones jack
5. AC adapter socket
6. MicroSD (TF card) slot
7. MicroHDMI out
8. Power on/off

**Caution!** Remember that mains voltage in the U.S. and Canada is 110V and 220-240V in most European countries. Please refer to the specifications table for the correct mains voltage and never attempt to plug a 110V device into 220V outlet and vice versa without using a converter.

## Installing a digital camera

Digital camera with LCD screen is installed on the third vertical tube of trinocular microscopes. Assemble and set up your microscope. Unscrew the upper part of the trinocular tube and then insert (screw in) the camera. Connect the supplied AC adapter to the camera and to the AC network. Press the power button and the camera will be ready to work.

## Embedded software

The embedded application is designed for using Android digital cameras with microscopes. With this application, you can change image parameters, measure objects as well as save images and videos for further use. In addition, you can perform particle analysis with its image processing feature. Like all Android applications, this application features a user-friendly multilanguage interface.

## Camera display interface

The screen is divided into two parts:

1. The left part is the **Preview Image**. It displays the current image provided by the camera.
2. The right part is the **Tool Panel**. It contains a number of tools for image capturing, changing parameters, measuring, etc. Use the right side of the screen to perform the following actions:

- Effect: Settings of the camera parameters
- Measure: Measuring tool
- Playback: Images and video playback
- Snap: Image capturing
- Record: Video recording
- Setting: Setting and changing parameters

## Capturing and playback tools

 Snap	Press the <b>Snap</b> button to take a photo. Select the image size in the <b>Settings</b> section.
 Record	Press the <b>Record</b> button to start video recording. To stop recording, press the same button once again. If the original size of the video is smaller than 1080p, then the size of the recorded video will be the same as the original. If the original size is bigger than 1080p, the recorded video size will be reduced to a size smaller than 1080p.
 Playback	Press the <b>Playback</b> button to open the last image captured or to play the last video recorded. The image will be opened in the <b>Image Analysis</b> window.

## Camera settings

The Effect panel allows you to edit the captured image, if necessary.

### Exposure

Auto Exposure: If you enable the checkbox next to “Auto Exposure”, the brightness of the image will be automatically adjusted.

Manual mode: You can set the exposure time and gain.

### Color temperature (White balance)

Auto mode: If you enable the checkbox next to “Auto WB”, the white balance will be automatically set.

Manual mode: For manual settings, use the “Color temperature” slider tool.

**Brightness**: Use the slider tool to adjust the brightness.

**Contrast**: Use the slider tool to adjust the contrast.

**Saturation**: Use the slider tool to adjust the saturation.

**Sharpness**: Use the slider tool to adjust the sharpness.



**Gamma**: Use the slider tool for the gamma correction.

**Flip image**: These icons are used to flip the image vertically or horizontally.

This icon is for saving the settings or resetting all of the settings to the default parameters. You can save up to four different parameter sets as well as return to the factory settings.

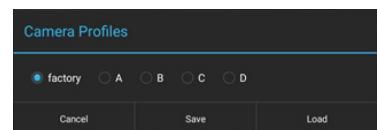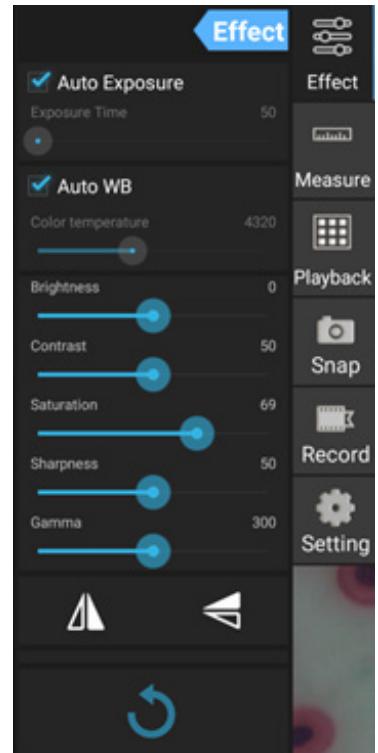
## Measurements

On the **Measure** panel, you can select a tool to conduct a measurement.


Before you start, you need to calibrate the ruler. You also need to calibrate it after changing the microscope objective and/or camera resolution.

### 1. Calibrations

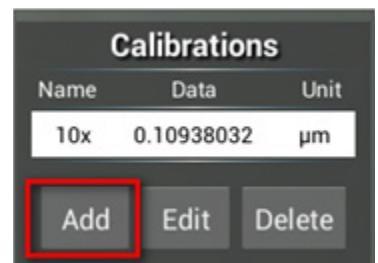Go to the **Measure** panel, and press **Add** button to start a new calibration process. Press the **Edit** button to make changes in previous calibrations.


Open calibration menu:

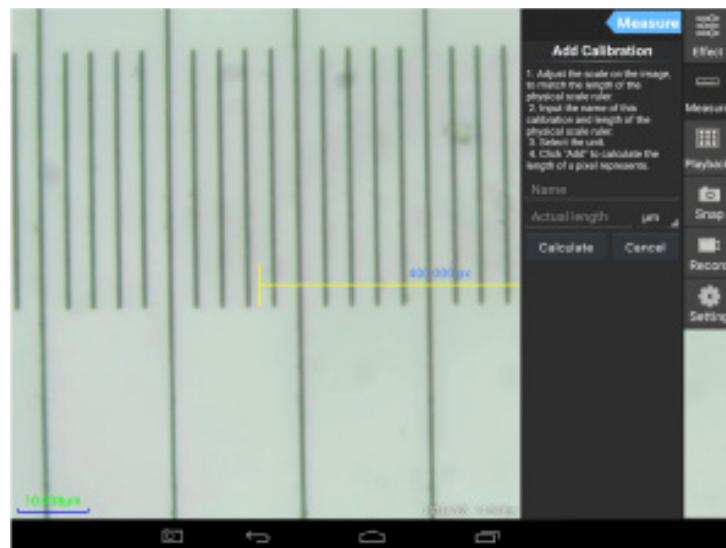Follow the instructions:

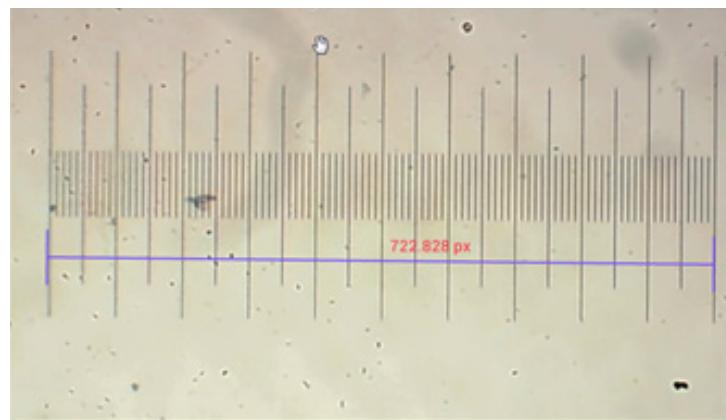Adjust the ruler to where its ends are on the physical ruler's scales. The ruler on the screenshot has 0.01mm scales, with each big grid equaling 10 $\mu$ m, and the whole ruler equals 1000 $\mu$ m.

Enter the name for the calibration and actual length of the ruler. In the example, the name of the calibration is "4x", which corresponds to the magnification value of the chosen eyepiece, and 1000 $\mu$ m is the actual length of the chosen ruler.

Press **Calculate** to get the calibration value for the current eyepiece and image size. Press **Add** to save the calibration.

The left screenshot shows the 'Add Calibration' dialog. It contains instructions, a magnification input field ('4x'), an actual length input field ('1000 μm'), and 'Calculate' and 'Cancel' buttons. The right screenshot shows a 'Calibrations' table with two entries:

Name	Data	Unit
4x	1.3842217	$\mu$ m
10x	0.10938032	$\mu$ m

Buttons for 'Add', 'Edit', and 'Delete' are located at the bottom of the calibration list.

## 2. Measurement tools

You can use a wide range of measurement tools (rulers).

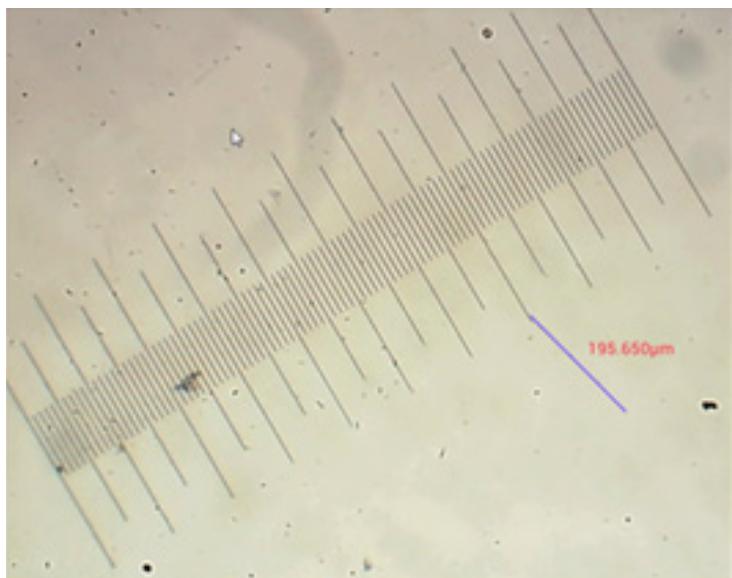
Icon	Measurement tool	Action
	Point counting	Set the starting point of measuring on the photo
	Line	Measure the distance between two points
	Rectangle	Measure the width, length, and area of the rectangle
	Circle	Measure the area of a circle
	Cross	Measure the cross hair
	Angle	Measure the angle
	Two circles	Measure the distance between the centers of two circles
	Perpendicular	Measure the length of a perpendicular line
	Concentric circles	Measure the radius of two concentric circles
	Text annotation	Add a comment to an image
	Option	Change the width or the color of the ruler's scale, and the size or the color of the text
	Export	Export the image with measure marking
	Delete	Delete the chosen ruler
	Delete	Delete all

## 3. Use of the measuring ruler

Choose a calibration (the chosen option will be displayed as a black print on the white background).


Calibrations		
Name	Data	Unit
4x	1.3842217	µm
10x	0.10938032	µm
<a href="#">Add</a>	<a href="#">Edit</a>	<a href="#">Delete</a>

Choose a ruler from the side menu, and it will appear in the image.

Use the measuring ruler to measure the physical ruler.

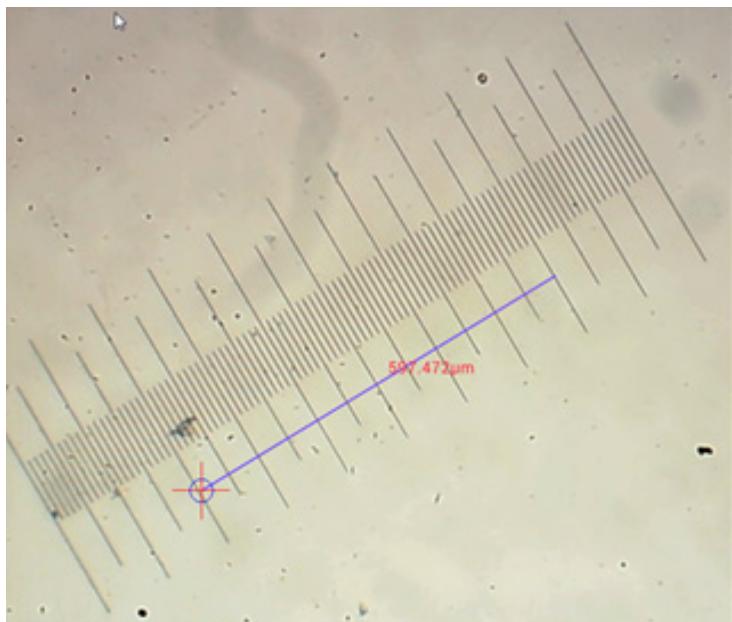
In this example, 6 large grids were measured: the measuring ruler shows  $597.472\mu\text{m}$ , which confirms the correctness of the measurements.

## Image analysis


The images are analyzed with the Android application YWImage, which uses OpenCV (Open Source Computer Vision Library). The application version 1.0 supports such functions as gray scale, contrast, binarization, and particle analysis.

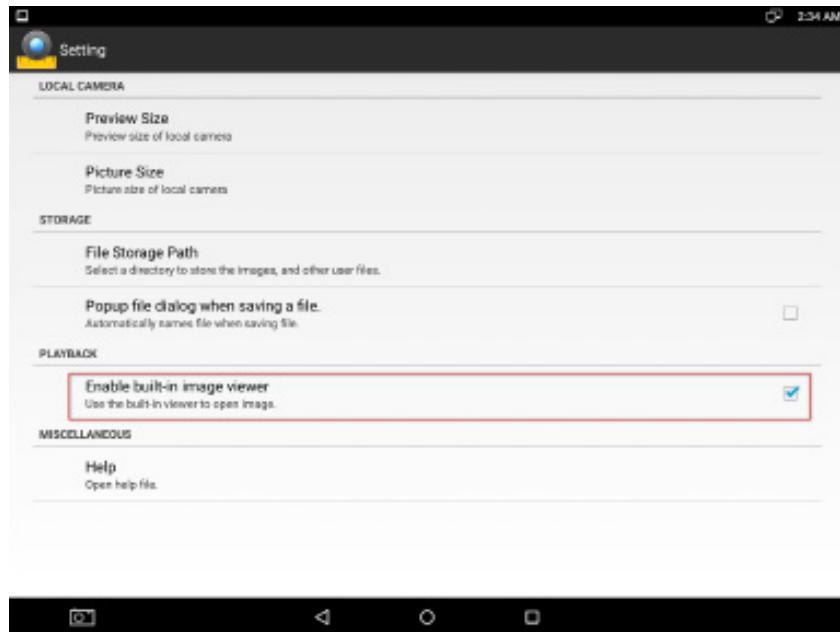To enable this feature, select the **Enable built-in image viewer** option in the settings dialog box.


Let us consider an example of particle analysis.

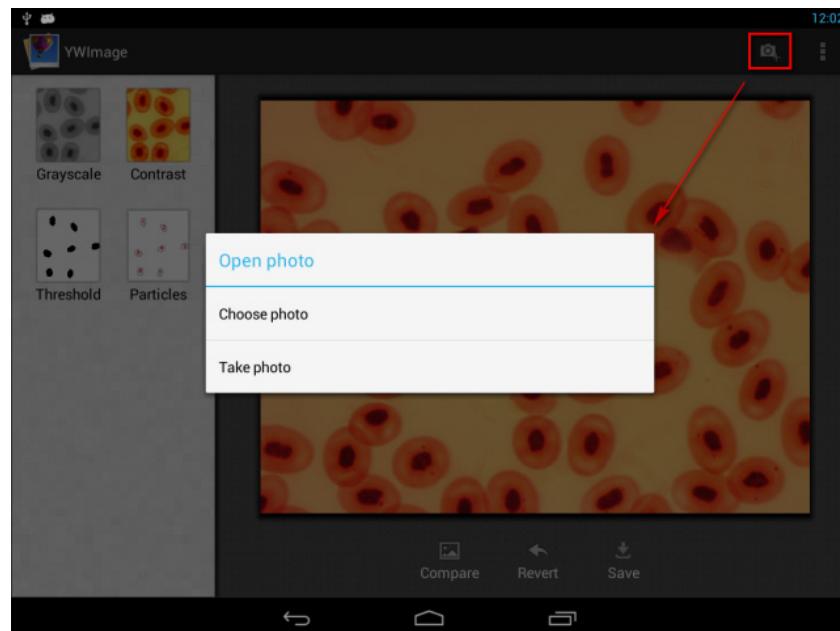
Image uploading. Click the camera icon in the upper right corner of the screen. You can select an existing image or take a picture with the camera.

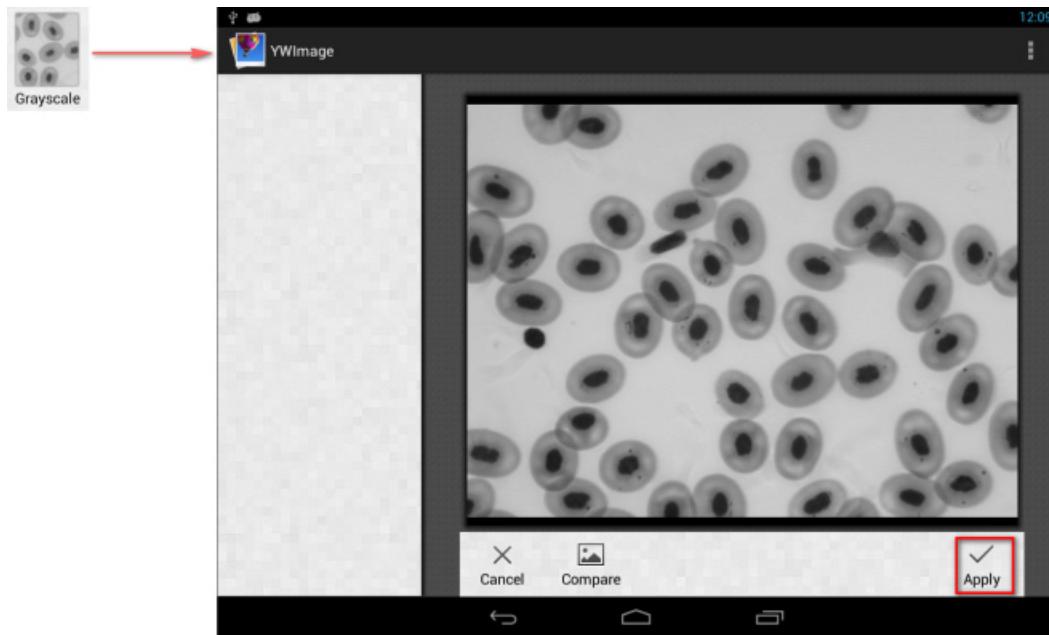Convert a color photo to Grayscale mode. Choose **Grayscale** from the left menu and press **Apply** to confirm converting.

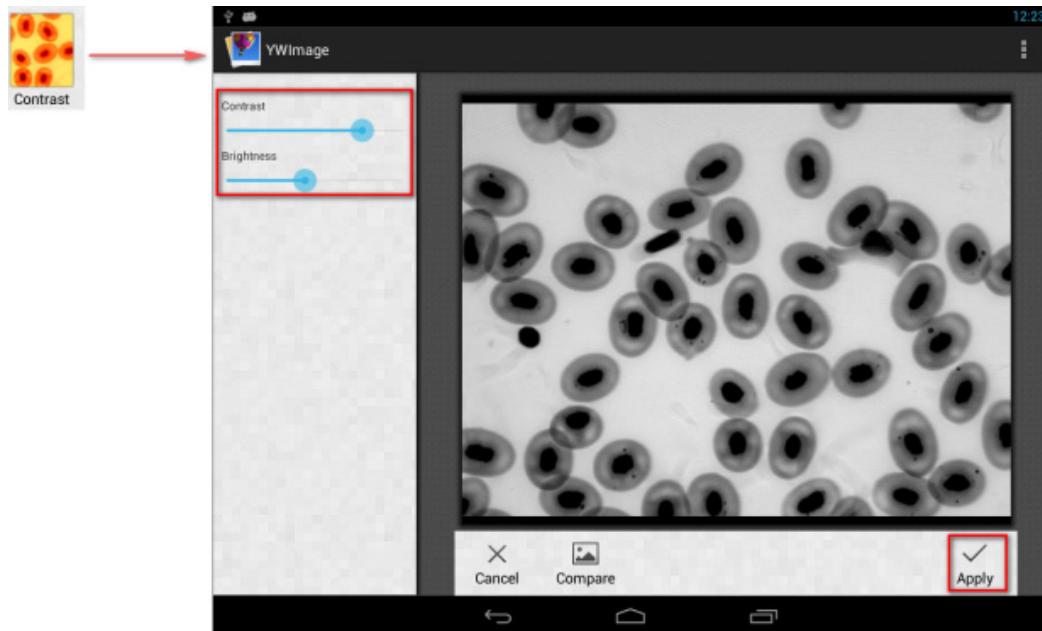
Contrast and brightness adjustment. Adjust the contrast and brightness of the black and white image, increasing the difference between objects and the background.


Binarization. Using the binarization function, you can convert a color image or an image in grayscale to two-color (black and white) by setting the minimum and maximum thresholds with which the brightness of each pixel will be compared.

All pixels within the threshold range will be highlighted in red. After executing the binarization, the RGB-value of the red areas will be (0, 0, 0), the RGB-value of the other sections will be (255, 255, 255).

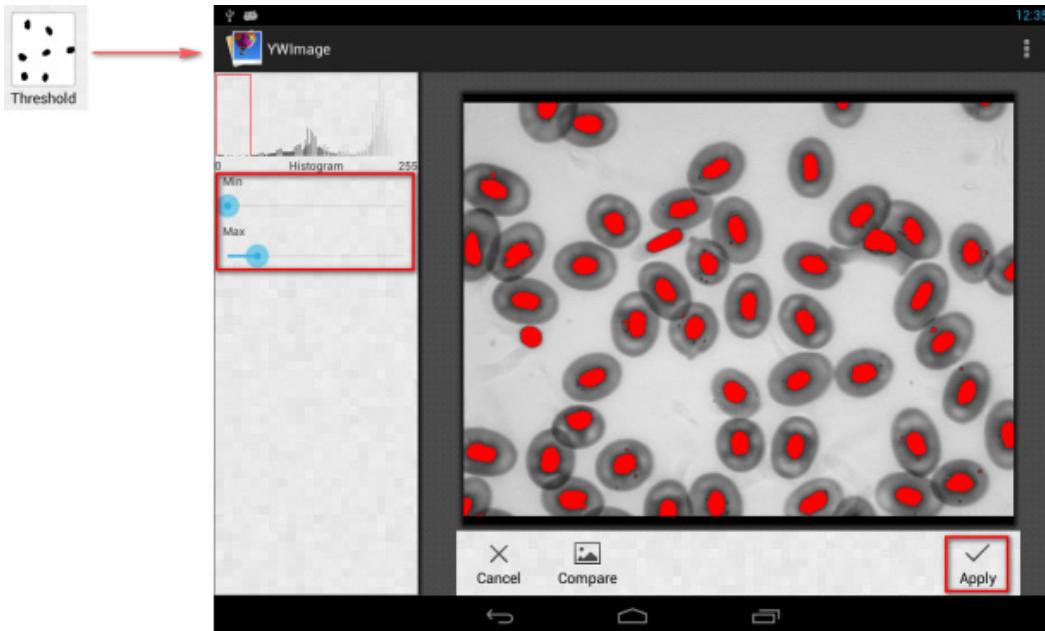**Note:** Red areas should not overlap.

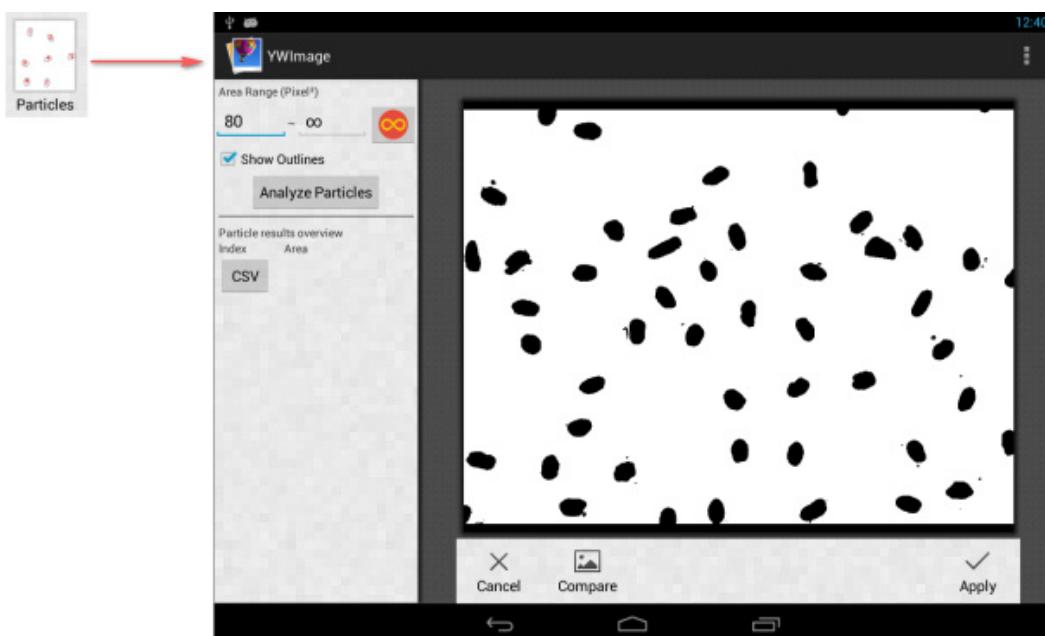Particle analysis. After executing the binarization, open the particle analysis tool.

First of all, if some noises were not deleted during binarization, you need to specify a particle size range. The designation of the particle size range will make the result more accurate.

Press the button , to set the maximum value to infinity. Select the Show Outlines check box to show the boundaries and number of particles after the analysis. Click **Analyze Particles** to begin the analysis. The result will be shown below.

Click the **CSV** button to export the data to the report file, which will be saved here by default:  
/mnt/sdcard/Image/Reports

## Specifications

LCD display	9.4" color touch screen
Megapixels	5
Maximum resolution (for still images)	2048x1536px
Sensor	1/2.5"
Pixel size	2.2x2.2µm
Frame rate	15 frames per second
Exposure control	auto/manual
Video recording	+
Image format	*.jpg
Video format	1080p, *.3gp
White balance	auto/manual
Output	USB 2.0 (2pcs.), microHDMI, Wi-Fi, TF card slot
Power supply	12V/2A; via AC adapter
Software features	measuring, brightness, exposure time, particle analysis, etc.
Mounting location	third eyepiece tube of a diameter of 23.2mm
OS	Android 5.1 (multilingual)
Embedded memory	4GB
Ability to connect additional equipment (not included in the kit)	support of microSD cards with capacity up to 32GB; monitor/TV (with a HDMI port); flash disk, mouse, keyboard (with a USB port); headphones (3.5mm)
The kit includes	digital camera with LCD screen, AC adapter

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The manufacturer reserves the right to make changes to the product range and specifications without prior notice.

## Care and maintenance

- Never look at the sources of bright light or lasers through the camera or the microscope: it will cause **PERMANENT DAMAGE TO YOUR EYES!**
- Do not disassemble the camera on your own.
- Keep the camera away from moisture; do not use it in rainy weather.
- Keep the camera away from shock or excessive pressure.
- Do not overtighten the locking screws.
- Keep the camera away from hazardous environment, home and car heaters, incandescent lamps or open fire.
- When cleaning the lenses, first blow any dust or debris off the surface or wipe them off with a soft brush. Then wipe the lens with a soft cleaning cloth slightly damp with spirits or ether.
- Seek medical advice immediately if a small part or a battery is swallowed.

## Levenhuk International Lifetime Warranty

All Levenhuk telescopes, microscopes, binoculars, and other optical products, except for their accessories, carry a **lifetime warranty** against defects in materials and workmanship. A lifetime warranty is a guarantee on the lifetime of the product on the market. All Levenhuk accessories are warranted to be free of defects in materials and workmanship for **six months** from the purchase date. The warranty entitles you to the free repair or replacement of the Levenhuk product in any country where a Levenhuk office is located if all the warranty conditions are met.

For further details, please visit: [levenhuk.com/warranty](http://levenhuk.com/warranty)

If warranty problems arise, or if you need assistance in using your product, contact the local Levenhuk branch.

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Two screenshots of the "Measure" dialog box from the software. The left screenshot shows the "Add Calibration" section with the following text:

1. Adjust the scale on the image, to match the length of the physical scale ruler.  
2. Input the name of this calibration and length of the physical scale ruler.  
3. Select the unit.  
4. Click "Add" to calculate the length of a pixel represents.

Below this, there is a text input field with "4x" selected as the unit. Underneath the input field, the value "1000" is entered into a dropdown menu next to the unit "μm". At the bottom are two buttons: "Calculate" and "Cancel". The right screenshot shows the "Calibrations" section of the dialog, displaying a table of calibration data:

Name	Data	Unit
4x	1.3842217	μm
10x	0.10938032	μm

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Two screenshots of a digital microscope software interface. The left screenshot shows the "Add Calibration" dialog box. It contains four steps: 1. Adjust the scale on the image, to match the length of the physical scale ruler. 2. Input the name of this calibration and length of the physical scale ruler. 3. Select the unit. 4. Click "Add" to calculate the length of a pixel represents. Below these steps are input fields: "Name" (text box with "4x"), "Actual length" (text box with "1000"), and "Unit" (dropdown menu set to "μm"). At the bottom are "Calculate" and "Cancel" buttons. The right screenshot shows a "Calibrations" table. It lists two entries: "4x" with a value of "1.3842217" and "Unit" "μm", and "10x" with a value of "0.10938032" and "Unit" "μm". At the bottom are "Add", "Edit", and "Delete" buttons.

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Two side-by-side dialog boxes from the software. The left dialog is titled "Add Calibration" and contains instructions for calibrating a 4x magnification image. It includes a text input field for "Name" containing "4x" and a dropdown menu for "Unit" set to "μm". The right dialog is titled "Calibrations" and lists two entries: "4x" with a value of "1.3842217" and "Unit" "μm", and "10x" with a value of "0.10938032" and "Unit" "μm". At the bottom of the right dialog are three buttons: "Add", "Edit", and "Delete".

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

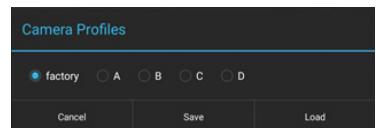
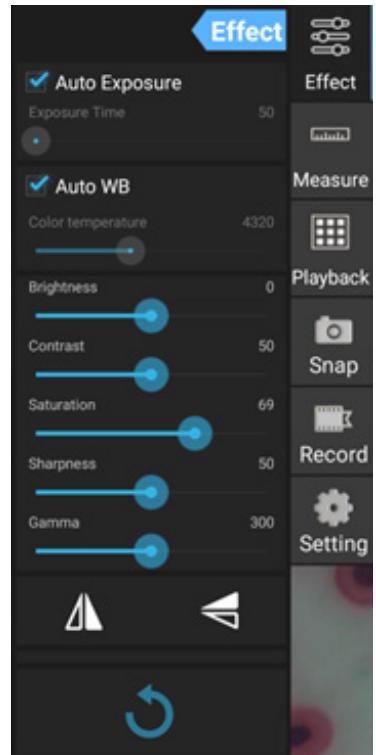
Calibrations		
Name	Data	Unit
10x	0.10938032	µm
<b>Add</b>	<b>Edit</b>	<b>Delete</b>

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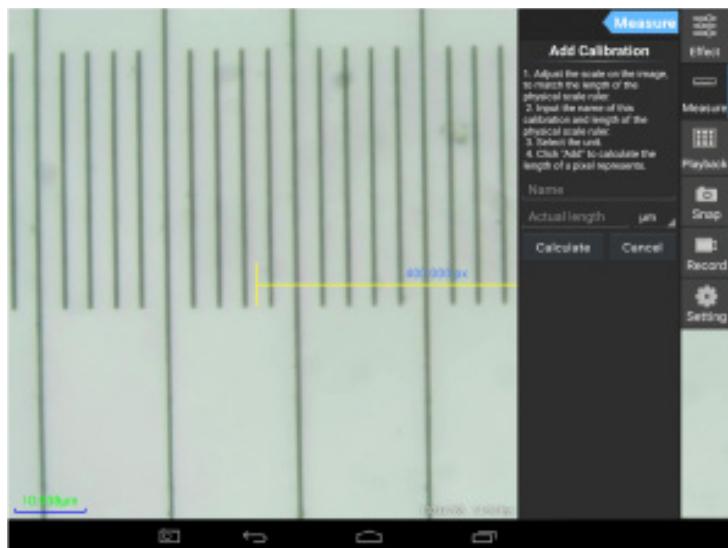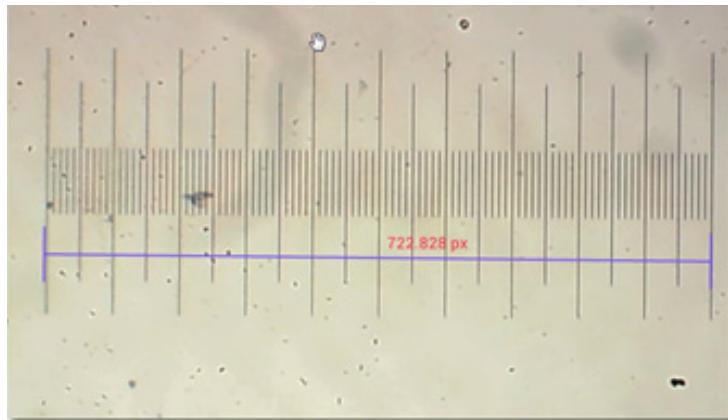Two screenshots of a digital microscope software interface. The left screenshot shows the "Add Calibration" dialog box. It contains instructions, a "Name" field with "4x", a "Data" field with "1000", a "Unit" dropdown set to "μm", and buttons for "Calculate" and "Cancel". The right screenshot shows the "Calibrations" table with two entries: "4x" and "1.3842217 μm" and "10x" and "0.10938032 μm". Below the table are buttons for "Add", "Edit", and "Delete".

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Two screenshots of a calibration dialog box. The left screenshot shows the "Add Calibration" steps: 1. Adjust the scale on the image, 2. Input the name of this calibration and length of the physical scale ruler, 3. Select the unit, 4. Click "Add" to calculate the length of a pixel represents. It also shows input fields for "Name" (4x), "Actual length" (1000), and "Unit" (μm). The right screenshot shows a "Calibrations" table with two entries:

Name	Data	Unit
4x	1.3842217	µm
10x	0.10938032	µm

With buttons for "Add", "Edit", and "Delete".

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Levenhuk® is registered trademark of Levenhuk, Inc.
© 2006-2025 Levenhuk, Inc. All rights reserved.
Levenhuk Inc. (USA): 928 E 124th Ave. Ste D, Tampa, FL 33612, USA,